



Ultra-sensitive Raman microscopy and endoscopy tools for pre- and intraoperative imaging


EPIC Meeting at NKI December 2019

Bodo Richter, APE GmbH Berlin





Raman Spectroscopy

- Chemical information is derived by probing vibrational states of molecules
- Finds wide applications, e.g. in material science
- Downside: rather small cross section => low intensities and sensitivity

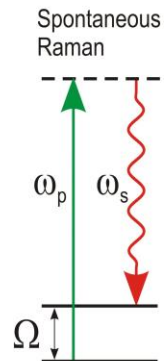


Stimulated Raman Spectroscopy/Microscopy

- Significant ($\times 10^6$) increase in sensitivity by using two laser beams
- Stimulated Raman Spectroscopy
- Vibrational excitation is drastically amplified when energy difference between Pump and Stokes matches the vibration of chemical bond of interest (e.g. CH_2)
- Images can be generated by scanning the lasers beams in a scanning microscope and detecting intensity changes at each pixel

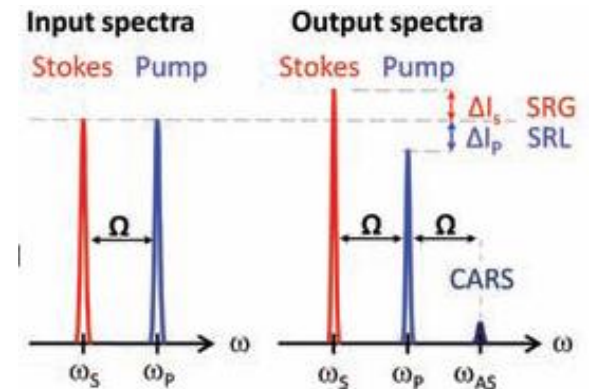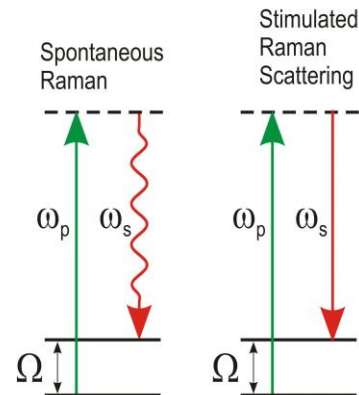

How can Stimulated Raman
Microscopy be used in
Histology?

Histology: H&E staining


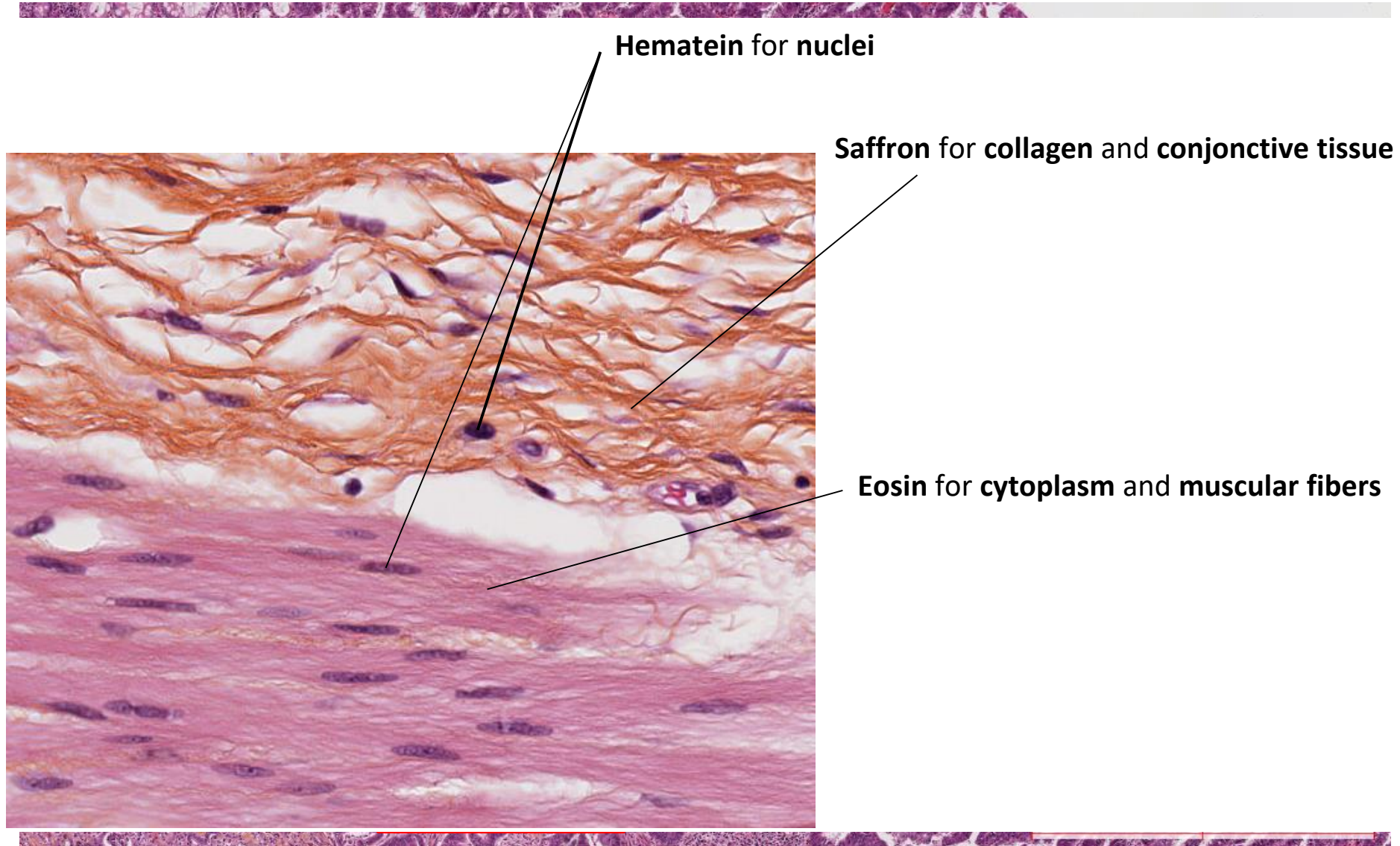


Histology: H&E staining

- Well established procedure
- Cutting – Fixing – Staining – Slicing
- Time consuming
- Could it be improved?

Requirement from the medical doctors - **Neurology**



- Intra-operative imaging of tumour tissue
'Cutting' has functional consequences
- Not necessarily H&E (Haematoxylin and Eosin) staining (5ALA)

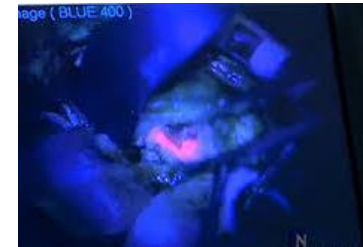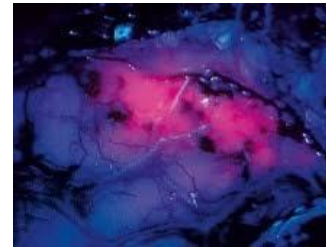

- Intra-operative means imaging **before** removing, ideally in a flexible probe

Requirement from the medical doctors - Neurology

- **BUT** a first step providing ex vivo H&E diagnostic is valuable
- Also neurobiologist say that the immediate diagnostic after brain biopsy is valuable

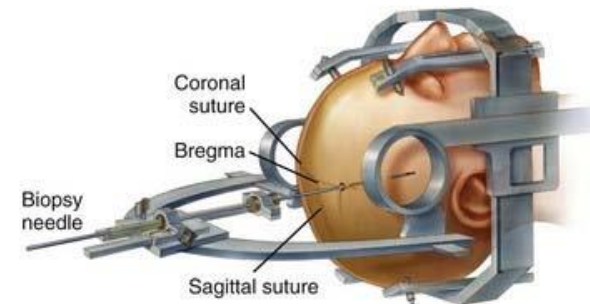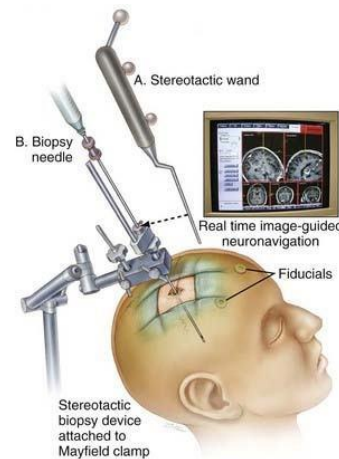

Requirement from the medical doctors – Gastric System

- Although gastric system removal is less critical than brain there are situations where instant histology would save lives
 - The detection of peritoneal metastasis during a gastric surgery
 - The clear identification of low and high grade epithelial dysplasia taking place at the gastroesophageal junction
- Flexible CRS probes can be inserted into the user port of commercial endoscopes



Goal




Is it possible to do:

- Label free histology with H&E quality?
- Instantaneous to be compatible with in vivo sample?
- Perform label free histology in an endoscope to access deep tissue and intra-operative diagnostic?

Stimulated Raman Histology (SRS)

Orringer et al., **Nat. Biomed. Eng.** **1**, 0027 (2017) – SRH in the brain

B. Sarri et al, **Scientific Reports** **9**, 10052 (2019) – SRH in the gastric system

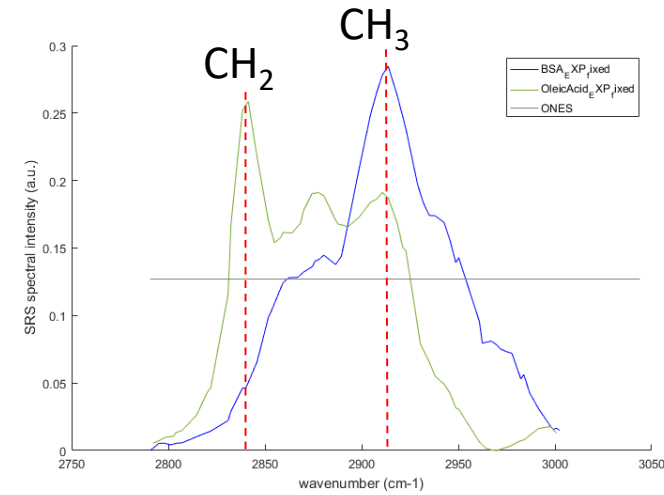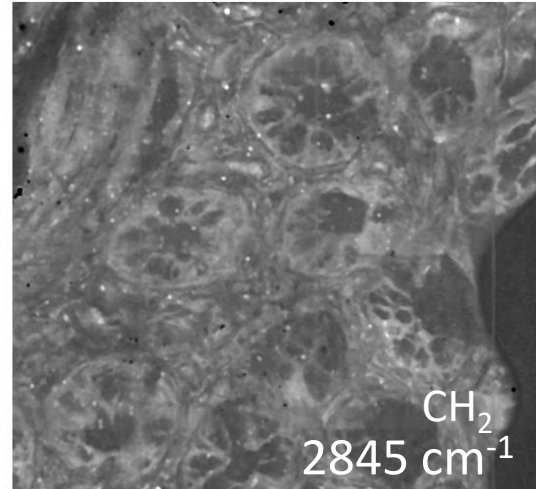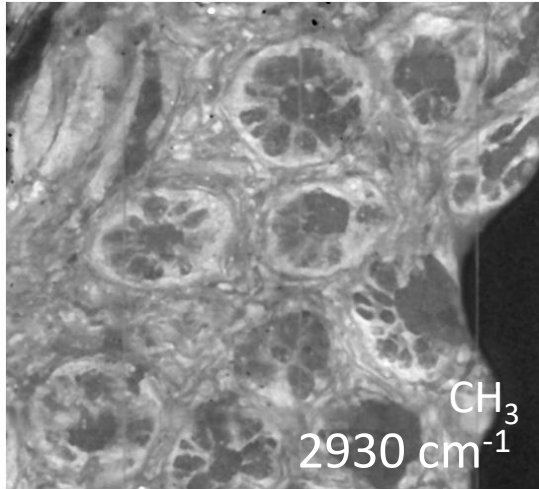


Fresh tissue


Diagnostic: SRH versus H&E :

Colon

Pancreas

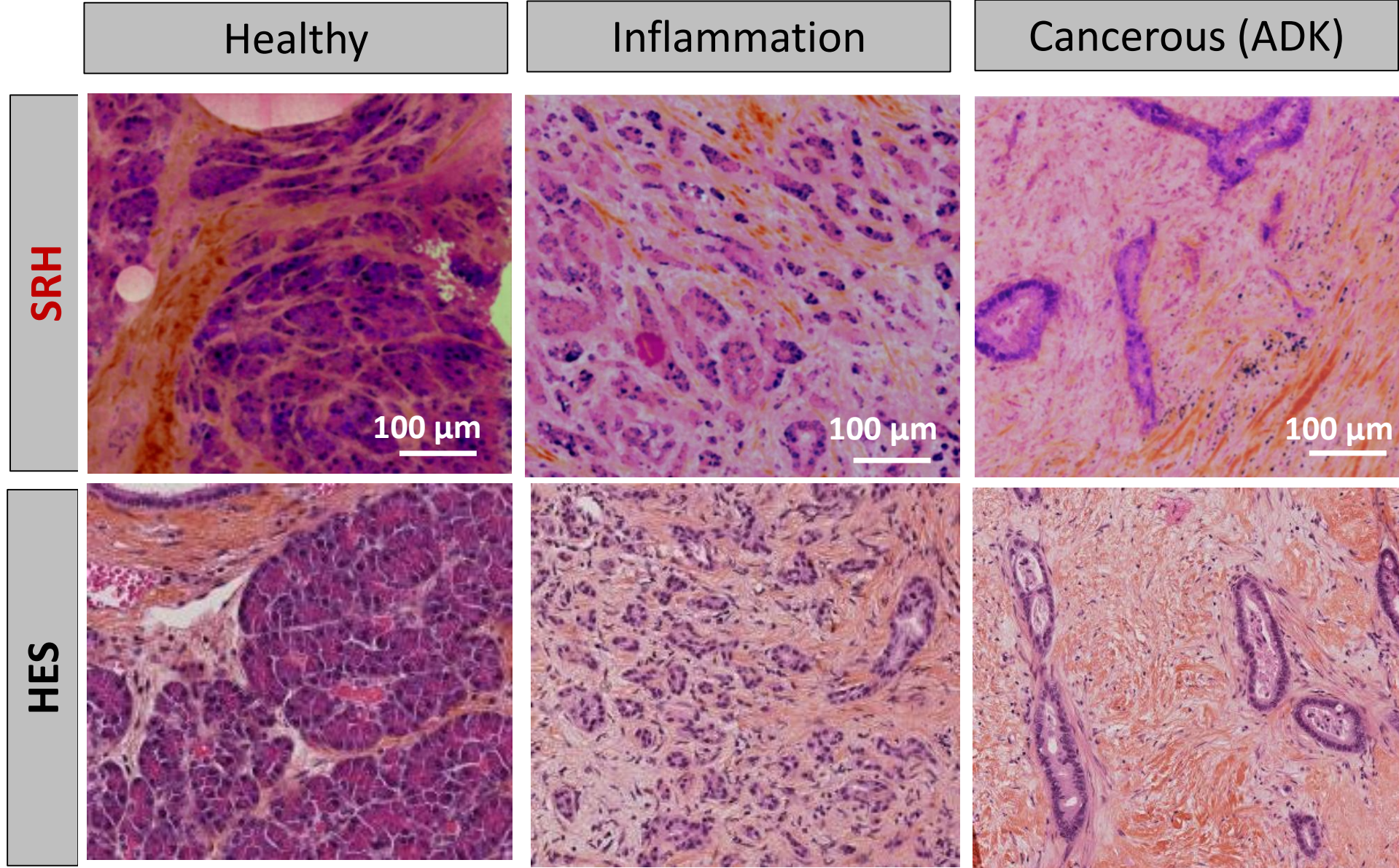

SRH Intraoperative context

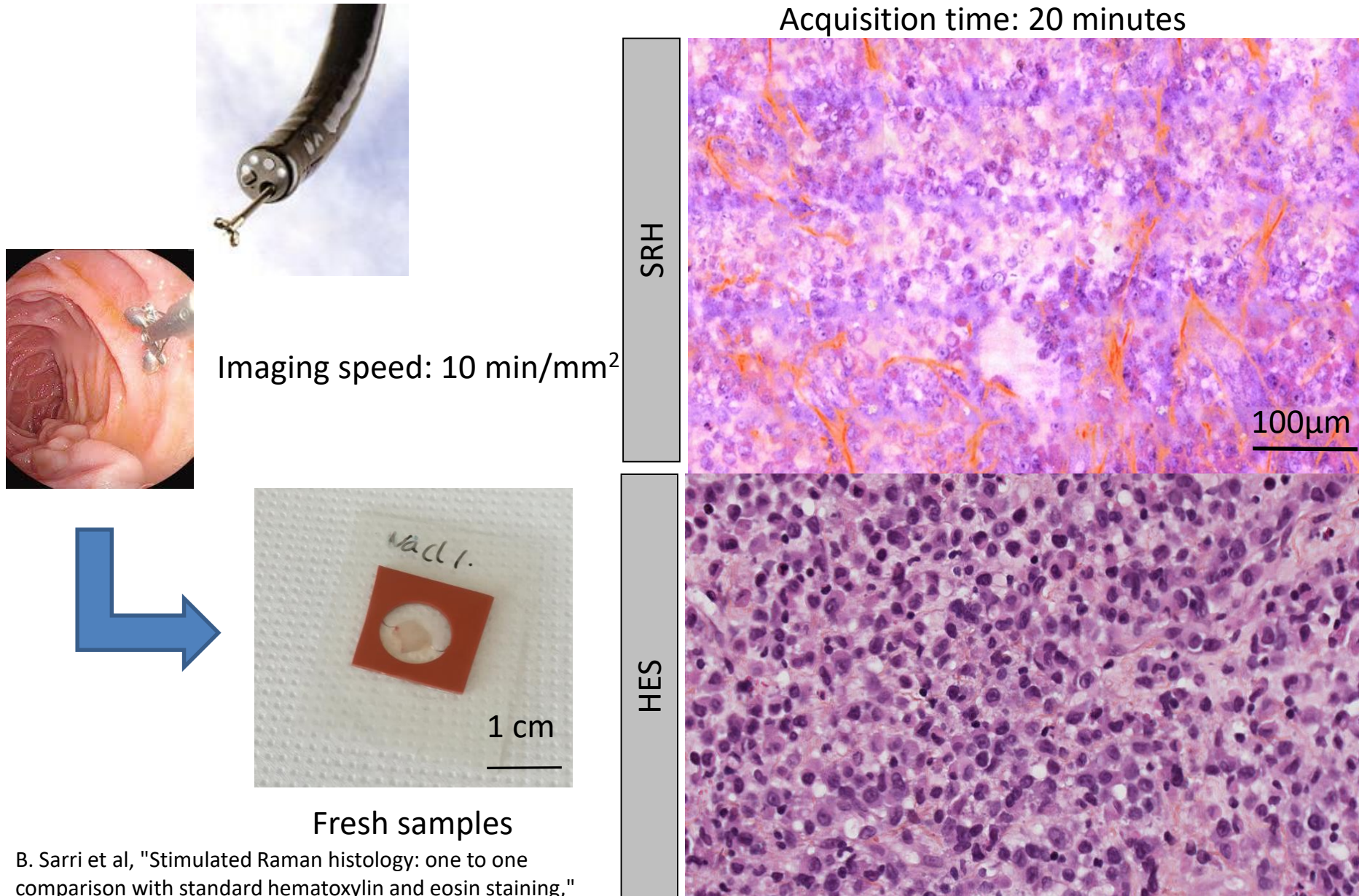

B. Sarri et al, "Stimulated Raman histology: one to one comparison with standard hematoxylin and eosin staining," Biomed. Opt. Express 10, 5378-5384 (2019).


Conclusion

- Stimulated Raman Histology (SRH) can provide almost equal information as H&E-staining
- Information is available much faster => important especially during operation
- SRH can provide additional information due to improved z-resolution; less distortions from cutting
- SRH can be applied to various types of human tissue
- Compatible with general workflow in operatory room
- Further clinical studies necessary to show benefit of technology

Outlook: Coherent Raman imaging system



- **Stimulated Raman histology**
- Coherent Raman imaging (CARS)
- 2-photon imaging
- Second harmonic imaging
- User friendly interface
- Millimeter field of view
- Sub-micron resolution
- 3D imaging capability
- 1mm² image in <60s

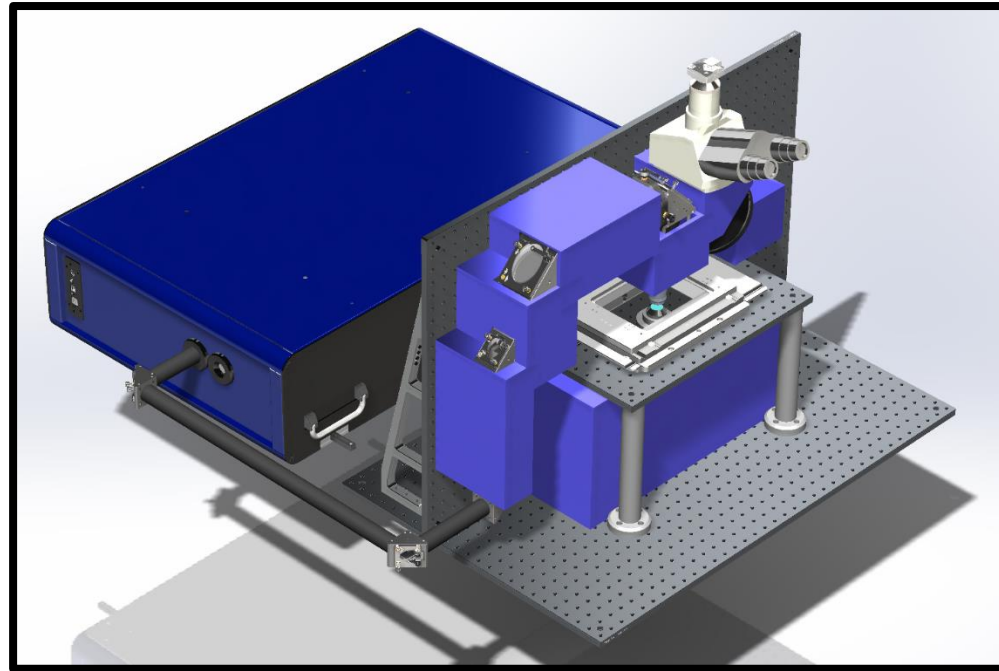

Available in 2020!

This presentation was presented at EPIC Meeting on Photonics for Cancer Diagnostics and Treatment 2019

HOSTED BY



SILVER SPONSORS



EU initiatives funded by
www.photonics21.org



PHOTONICS PUBLIC PRIVATE PARTNERSHIP

BRONZE SPONSORS

